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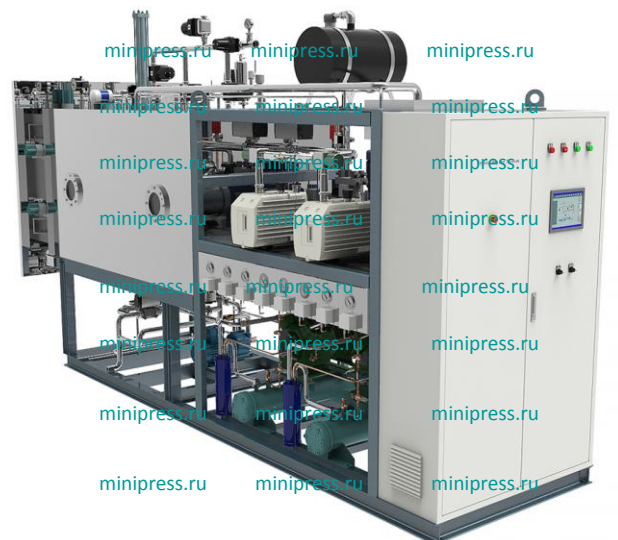
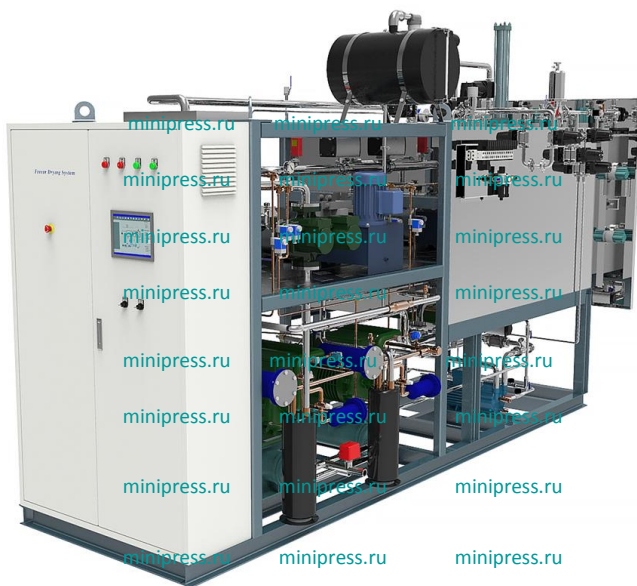
FABRICANTE - CHINA

SPECIFICATIONS

We represent the largest Chinese manufacturer specializing in sublimation equipment in Russia and Europe.

By order of the customer, freeze drying can be designed and manufactured with any number of shelves and usable area.

 www.minipress.ru

ADDITIONAL INFORMATION

The process of operation of this equipment consists in placing in the freeze drying chamber vials of liquid material, followed by obtaining dried powder. Pre-rinsed and dried glass penicillin bottles are used in which the material is poured. Then the vials are sealed with special rubber stoppers having side slits, in the part that enters the neck of the vial. Capping is not done until the end, but by 50%, only to fix the rubber stopper in the neck of the bottle. Subsequently, through the slotted vacuum, the liquid is pumped out of the vial. In the final stage, after complete dehydration of the material, the cork is completely capped by hydraulic moving shelves. Thereby, the drying takes place in a vacuum medium, and then the material is finally packed in a vacuum. This gives (in contrast to other types of such drying) to keep the material in a vacuum for 1-2 years. After the vials are evacuated, the chamber can be opened and an automatic or semiautomatic fixation of the rubber stopper with an aluminum cap can be made. Caps are put on the neck of the bottle and sealed. According to the manufacturer of the equipment, with this technology it is possible to completely restore the life of the material after dehydration at the opening of the vacuum package within 4-5 hours. Storage in a vacuum of the material is possible at a temperature not exceeding +36 degrees C. Since in the process of dehydration (removal of water) the equipment dehydrates the material, then the recovery occurs with distilled water. The water to be removed is frozen in the chamber, then drained after defrosting and it is possible to control how many ml left from the dried product in order to subsequently restore just such a quantity of water. This is an important point. According to the manufacturer, not all materials can be turned into powder, some biological materials become dehydrated into a jelly-like substance, which is also reconstituted with distilled water. The equipment manufacturer has a large laboratory that, together with the pilot production, can conduct any research of your material, obtain samples of the products in dried form and packed in vacuum for further research by customers.

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